

## **Hart's Stain of Paraffin Sections**

1. Americlear for 15 min.
2. Deparaffinize - 100%, 95%, 80%, 70%, 50%, 30% ETOH, H<sub>2</sub>O (10 dips each)
3. Stain in Rescorcin Fuchsin (Hart stain) for 1-2 hours (can be reused)
4. Rinse with tap water 5 min.
5. Dehydrate - 30% ETOH, 50%, 70%, 80%, 95%, 100%, 100% (10 dips each)
6. Americlear for 5 min.
7. Coverslip with Permount

### **Rescorcin Fuchsin Working Soln.**

- 10 ml Rescorcin Fuchsin stock (*can be purchased from Polyscientific.*)
- 100 ml 70% ETOH
- 2 ml Conc. HCl